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Thành phần:

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Product: 5PRP Kit V1

Components:

Spinal anesthesia needle; 5 mL Luer-lock Syringe; Filter; Connector; 10 mL Luer-lock Syringe; Alcohol pad; Kim Vacuum blood collection needle; Needle holder; Bandage; PRP tube; The PRP storage tube.

Reference number: 376.

Amount: Kit/box

Purpose of use:

For research or manufacturing use only.

5PRP Kit V1 Kit is used to isolate and activate platelet-rich plasma from 12ml of peripheral blood, umbilical cord blood, etc.

Summary of salient features

5PRP Kit V1 is a Kit for blood collection and manufacturing of activated platelet-rich plasma. This Kit draws blood using vacuum tubes and separates platelets by isopyclic centrifugation. PRP is activated through a filtration membrane, resulting in a final activated PRP product that does not contain any blood cells and has a high purity level. The product manufacturing process is done quickly with 1 single centrifugal step. 5PRP Kit V1 was researched and developed by Stem Cell Institute, HCMUS, VNUHCM, Viet Nam. 5PRP Kit V1 guarantees to be sterile (0 CFU) with endotoxin level ≤ 1 EU/mL.

Reviewed/tested Applications:

5PRP Kit V1 provides a fast and convenient method to collect blood and manufacture platelet-rich plasma from peripheral blood or umbilical cord blood, etc.

Storage conditions and shelf life:

Storage and transportation temperature: Room temperature.

Expiry date: following the nearest expiry date of the components

Procedure:

5. Make sure the product is intact and within the expiry date.
6. **Blood collection**
 - 6.1. Attach the Vacuum blood collection needle (butterfly needle) to the Needle holder.
 - 6.2. Clean the puncture site with an alcohol pad, then perform venipuncture.
 - 6.3. Advance the PRP tube into the Needle holder; at this point, blood will flow automatically into the tube.
 - 6.4. Collect blood until the tube is full (tube contains 10-12 mL of blood). Gently invert tubes to mix blood with an additive.
7. **PRP Production:**
 - 7.1. Place the PRP tubes containing blood into the centrifuge (place 1 more counterweight tube in a symmetrical position).
 - 7.2. Centrifuge the tubes at **1500g for 8 minutes**.
 - 7.3. Gently take out the PRP tubes from the centrifuge.
 - 7.4. Attach the 10 mL Luer-lock Syringe to a Spinal anesthesia needle. Remove supernatant on the top and leave **1,0-1,5 mL of plasma** at the

Regenedlab

Cellatist

Stem food
Your Genome And

Dermaloka

bottom (containing platelet pellet). *Please see the FactSheet to know about number of platelets in the tube.*

7.5. Suspend the platelet pellet in the plasma, and aspirate 3-5 times. This mixture is platelet-rich plasma (PRP).

8. Activation and sterilization of PRP:

8.1. Collect inactivated PRP into the 10 mL Luer-lock Syringe.

8.2. Attach the 10 mL Luer-lock Syringe to the Filter.

8.3. Attach the Connector to the other end of the Filter.

8.4. Attach the 5 mL Luer-lock Syringe to the other end of the Connector.

8.5. Strongly push the piston of the 10 mL Luer-lock Syringe so that inactivated PRP goes through the Filter and into the 5 mL Luer-lock Syringe (The product in 5 mL Luer-lock Syringe is activated PRP).

Precautions:

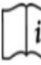
Do not use the Kit if the packaging is damaged or cracked, or if one of the Kit components is abnormal.

Troubleshooting:

Not applicable

Explanation of symbols and warnings

The symbols present on the Kit are explained below

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To purchase other products, please visit: <http://biomedmart.org>

For further information, please contact: contact@sci.edu.vn; sales@sci.edu.vn;

Regenedlab

Cellatist