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SẢN PHẨM: SPRP

Thành phần:

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Product: SPRP

Components:

Spinal anesthesia needle; 5 mL Luer-lock Syringe; Filter; Connector; 10 mL Luer-lock Syringe; Alcohol pad; Vacuum blood collection needle; Needle holder; Vacuum blood collection tube; Bandage; Plasma and platelets containing tube

Reference number: 170 (SPRP Kit); 190 (SPRP Kit V1)

Amount: Kit

Purpose of use:

For research or manufacturing use only.

SPRP is a kit used to isolate and activate platelet-rich plasma from peripheral blood, umbilical cord blood...

Description:

SPRP is the kit for blood collection and manufacturing of activated platelet-rich plasma. This kit draws blood using vacuum tubes and separates platelets using the centrifuge method. In particular, PRP is activated according to the mechanical principle, therefore, the final activated PRP product does not contain any blood cells and is highly purified. The SPRP kit was researched and developed by Stem Cell Institute, HCMUS, VNUHCM. . SPRP kit is tested as sterile (0 CFU) and has low endotoxin level (≤ 1 EU/mL).

Applications:

SPRP can easily and successfully collect blood and manufacture platelet-rich plasma from peripheral blood or umbilical cord blood...

Storage conditions and shelf life:

Storage and transportation temperature: Room temperature.

Expiry date: following the nearest expiry date of the components

Protocol:

Make sure the product is intact and within the expiry date.

5. Blood collection

- 5.1. Attach Vacuum blood collection needle (butterfly needle) to the Needle holder.
- 5.2. Clean the puncture site with an alcohol pad, then perform venipuncture.
- 5.3. Advance the Vacuum blood collection tube into the Needle holder. The blood will automatically draw into the tube.
- 5.4. Collect blood until tube(s) is (are) full (each tube contains 10-12mL of blood). Gently invert tubes to mix blood with an additive.

6. Centrifugation

- 6.1. Place Vacuum blood collection tube(s) into the centrifuge (place 2 tubes in a symmetrical position)
- 6.2. Centrifuge the tubes at **600g for 5 minutes** (if the centrifuge has brake option, turn it off before centrifuge the tubes)
- 6.3. Gently take out 2 Vacuum blood collection tubes from the centrifuge, and place on a 15 mL tube rack (the blood will be separated into 2 parts: blood cells are located at the

Regenmedlab

Cellatist

Stemfood

bottom and the yellow plasma layer containing platelets is on the top)

- 6.4. Use Spinal anesthesia needle to collect the platelet layer and plasma into the Tube containing plasma and platelets.
- 6.5. Place the Tube containing plasma and platelets in the centrifuge, and continue centrifugating at **900g for 15 minutes**.
- 6.6. After centrifugation, platelets settle at the bottom of the tube
- 6.7. Attach the 10 mL Luer-lock Syringe to a Spinal anesthesia needle. Remove the supernatant on the top and leave 1,0-1,5 mL of plasma at the bottom (containing platelet pellet).

Please read more information about concentration and dose of platelet harvested per tube in FactSheet

- 6.8. Suspend the platelet pellet in the plasma, and aspirate 3-5 times. This mixture is inactivated platelet-rich plasma (PRP)

7. Activation and sterilization of PRP:

- 7.1. Collect inactivated PRP into the 10 mL Luer-lock Syringe.
- 7.2. Attach the 10 mL Luer-lock Syringe to the Filter.
- 7.3. Attach the Connector to the other end of the Filter.
- 7.4. Attach the 5 mL Luer-lock Syringe to the other end of the Connector.

- 7.5. Strongly push the piston of the 10 mL Luer-lock Syringe so that inactivated PRP is filtered into the 5 mL Luer-lock Syringe (The product in 5 mL Luer-lock Syringe is activated PRP)

Cautions:

Do not use the Kit if the packaging is damaged or cracked, or if one of the Kit components is abnormal.

Troubleshooting:

Not applicable

Explanation of symbols and warnings

The symbols present on the Kit are explained below

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The symbols present on the product are explained below

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To purchase other products, please visit: <http://biomedmart.org>

For further information, please contact: contact@sci.edu.vn; sales@sci.edu.vn; kinhdoanh@sci.edu.vn